

Assembling guide for Re-Useable Jacket model BioStat



Overview of all components



1 step; put the SUB or SUF in the RUJ

Make sure it's centered



2 step; Place the Top ring with the flat side up



3 step; screw 3 wing nuts next to each other



4 step; screw the two support rod and the holder for the sky support, tightly. Be sure to retighten all six, to ensure it is closed



5 step; place the motor adapter




6 step; Place the nylon sprocket



7 step; Place the rubber part




8 step; Place the motor. Make sure to tighten the 4 bolts

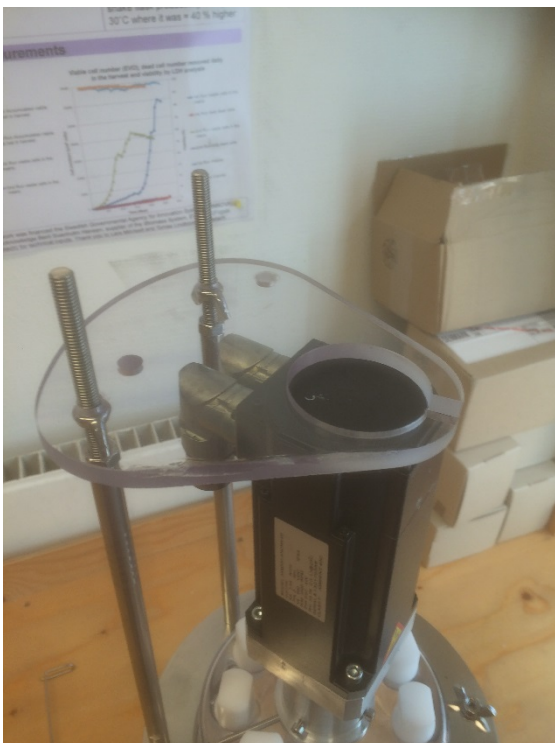

9 step; Place the motor holder



10 step; place the small round double sticky label on the motor

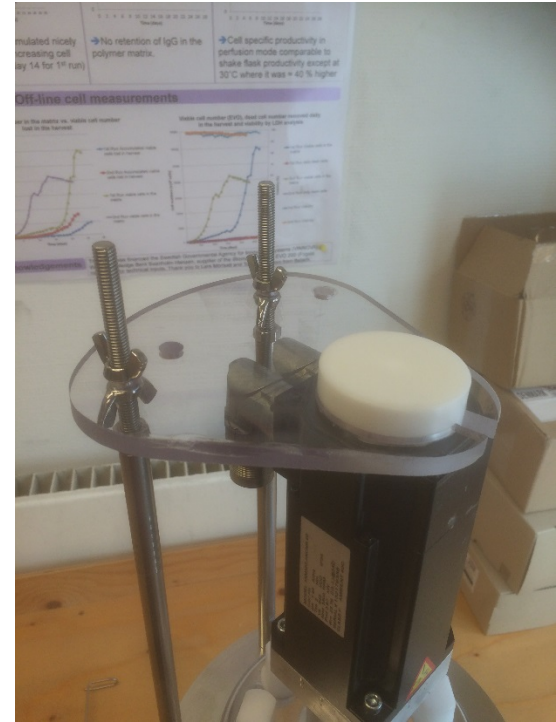

10 step; place the white nylon disk on the sticky label, in the center of the motor holder.



11 step; The RUJ is now ready for use