

# Assembly guide for RUJ/MST

This guide applies to the following number: 2286; 2287-xxx and 2288-xxx



## Item list



Number	Description
1	Main RUJ/MST
2	Your own SUB
3	Your own MST (not included) see: <a href="http://cercell.com/products/accessories/drive-system-s/magnetic-stirrer-table/">http://cercell.com/products/accessories/drive-system-s/magnetic-stirrer-table/</a>
4	4 pcs. M4 x 50 mm bolt
5	4 pcs. plastic dish with ID4,5
6	Your own adapter (not included) see: <a href="http://cercell.com/products/accessories/thermal-system-s/water-heat-cool-jacket-mst/">http://cercell.com/products/accessories/thermal-system-s/water-heat-cool-jacket-mst/</a>
7	RUJ cover ring
8	6 pcs. M6 wing nut
9	6 pcs. plastic dish with ID6,5



Start by unscrew the four bolts holding the support plate, and remove it from the MST.



Until the MST look like this



Place your reactor on the MST and align the holes in the MST with the ones in the RUJ



Place the ID4.5 dishes and screw the M4 x 50 on the RUJ. Make sure they're tighten firmly.





Make sure it tighten firmly



You're now ready to mount our own adapter



Make sure it tighten firmly



Careful you place SUB down in RUJ



Make sure it is center in the RUJ



Place the cover ring on the RUJ



Place the ID6.5 dish's and screw the wing nuts on the RUJ. Make sure they're tighten firmly.



You're now ready to use the RUJ, last step is to connection your own water hoses and motor